

Crosslinking

MASCOT



Mascot Server 2.7 and later can identify crosslinked peptides from MS/MS data. I'll show you how to use the new feature and discuss validating crosslinked search results.

Crosslinking experiments

- **Naturally occurring**
 - E.g. disulfide bond
- **Artificially created for elucidating:**
 - Protein tertiary structure
 - Protein complex topology
 - Conformational changes
 - Protein-protein interactions

Crosslinks can be roughly divided into naturally occurring and artificially created. An example of a naturally occurring crosslink is the disulfide bond. In most experiments, disulfide bonds are of no interest, so they are cleaved with a reducing agent prior to analysis. But, if the goal is to study protein folding, identifying the sites linked by a disulfide bond is crucial.

Crosslinks can also be created artificially by a chemical reaction. A crosslink can only be formed between sites that are adjacent in three-dimensional space. The maximum distance depends on the spacer arm length of the linker. So, the locations of the identified links can be used for constraining the three-dimensional shape of the molecule, which has obvious applications in elucidating protein tertiary structure, protein complex topology, conformational changes and protein-protein interactions.

Crosslinking experiments

- **Enzymatic digestion produces:**
 - Linear peptides
 - Crosslinked peptides: *alpha* linked with *beta*
- **Cleavable crosslink**
 - Cleaved prior to analysis, e.g. irradiation or CID
 - Leaves behind a fragment at linked site
- **Intact crosslink**
 - Survives MS/MS analysis
 - Spectrum has fragments from both peptides

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science

 **MATRIX
SCIENCE**

Whether natural or artificial, a typical bottom-up crosslinking experiment starts by digesting the crosslinked proteins. This produces two types of precursor: linear peptides and crosslinked peptides. A linear peptide either doesn't span a linkable site or if it does, the linker did not form a crosslink. Otherwise, the precursor is a crosslinked peptide, where an *alpha* peptide is linked with a *beta* peptide. The *alpha* peptide is conventionally the longer or more massive of the two.

How crosslinked peptides appear to a search engine like Mascot depends on the type of the linker.

Cleavable crosslinks are designed to be cleaved prior to analysis by various means like irradiation or CID. The link cleaves asymmetrically, so that the *alpha* and *beta* peptides carry linker fragments with different mass. The fragment is modelled exactly like a variable modification on a linear peptide. The linear peptide matches are then paired up by a postprocessing algorithm, which figures out which peptides come from the same crosslinked precursor.

Conversely, intact crosslinks survive MS/MS analysis. There is additional complexity to the crosslinked spectrum: *alpha* fragments that contain the linked site have a mass offset equal to the *beta* peptide mass plus the linker molecule. Vice versa for *beta* fragments. The search engine must be able to fragment crosslinked peptides *in silico*. However, the advantage is, no postprocessing step is needed.

The boundary between cleavable and intact is a bit fuzzy. If the cleaving efficiency

isn't 100%, there will be some precursors with an intact link. Similarly, an intact linker can fail to form a link, which leaves behind a linker fragment.

Introduced in Mascot Server 2.7

- **Model the experiment with a *crosslinking method***
 - Linker, linkable sites, directionality, crosslinked products (intralink, interlink, looplink, monolink), protein accessions, filters
- **Plus standard search parameters**
 - Applies to both linear and crosslinked peptides
- **Mascot ships with 14 linker definitions (Unimod), or create your own**

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science



Because of the great variety and complexity in crosslinking experiments, we've chosen an approach similar to quantitation. We want to keep the software universal and configurable, and not tied to a specific experimental workflow.

The crosslinking experiment is modelled with a crosslinking method. The method defines which linker is used, which linkable sites are considered, linker directionality, what are the expected crosslinked products and which protein or proteins are being linked. There are also a few filters available.

There are four possible crosslinked products. An intralink is an intact crosslinked peptide within the same protein, while an interlink is an intact link between two different proteins. A looplink is an intact link within a peptide. Finally, monolinks model cleavable linker fragments as well as dead-end or quenched links where only one end is attached to a peptide. I'll show examples of crosslinked products later.

The standard search parameters like enzyme, fixed mods and variable mods are applied to both linear peptides and crosslinked peptides.

The linker definitions are standard Unimod entries. Mascot currently ships with 14 definitions: 8 for intact crosslinks and 6 for cleavable crosslinks. You can easily create your own definition for any linker chemistry.

Data preparation - Peak picking

- **Mascot Server only searches for 1+ and 2+ fragment ion series**

- De-charge the fragment ion peaks to MH+
- Or supply the peak charge states
- For example, if you have a Thermo instrument use Mascot Distiller and the `prof_prof.ThermoXcalibur.opt` processing options

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science




There are several important factors to take into account when carrying out peak picking for an intact crosslinked dataset. You'll typically be dealing with higher charge state precursors to handle the larger masses of the linked peptides, and the MS/MS spectra are inherently complex, chimeric spectra with fragments from the alpha and beta peptides.

Mascot Server only searches for 1+ and 2+ fragment ion series. Crosslinked peptides are likely to have higher charge state precursors and many of the fragment ions generated will similarly have higher charge states so can't be matched (or worse, might result in false positive matches). For crosslinked datasets, you should either de-charge the fragment ion peaks to MH+ or supply the fragment peak charge states in the peak list as the third column.

Mascot Distiller can do this, but if your MS/MS scans are saved as centroids, rather than profile, then you need to use the uncentroiding options. For example, if you have a Thermo instrument and use `prof_prof.ThermoXcalibur.opt` processing options Distiller will uncentroid the data before processing. Peak picking time will be longer than using the default `ThermoXcalibur.opt` but the search results will be much improved.

Data preparation - Peak picking

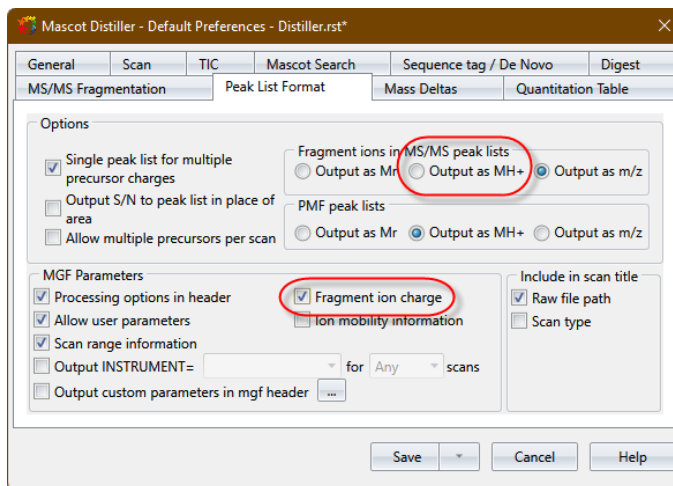

MASCOT

: Crosslinking

© 2020-2023 Matrix Science

**MATRIX
SCIENCE**

If using Mascot Distiller don't forget to set the peak list format to save the MS/MS peak lists output as MH+ or include the Fragment ion charge. If you make this change while a file is open it will only affect the current file. If you make the change while no files are open it will become the default both for when Distiller is used directly and for peak lists produced by Mascot Daemon using the Distiller data import filter.

Interlinked HOP2-MND1 complex

- Rampler et al: “Comprehensive Cross-Linking Mass Spectrometry Reveals Parallel Orientation and Flexible Conformations of Plant HOP2-MND1”, J Prot Res 14(12):5048-5062, 2015
- PRIDE: PXD001538
- Purified *A. thaliana* HOP2-MND1 complex
- DSS/BS3 linker, biological replicate 1

I'll illustrate intact crosslinking with this data set from the Mechtler lab. The data is available as the PRIDE project PXD001538. The study uses a purified HOP2-MND1 complex from *Arabidopsis thaliana*, which is crosslinked using three different linkers. I've chosen the biological replicate 1 linked with disuccinimidyl suberate (DSS). DSS has two identical reactive groups, so it's homobifunctional. It's amine reactive and creates intact crosslinks between lysines and protein N-term. The study was designed so that there are only lysine-lysine links.

Interlinked HOP2-MND1 complex

- **Peak pick in Distiller**

- Crosslinked precursor charge often 3+ or more, so output fragment masses as MH+

- **Set up a crosslinking method**

- Linker specificity: Xlink:DSS (K)
- Monolinks: enable W (156 Da)
- Accessions: HOP2_ARATH, MND1_ARATH
- Scope: enable InterLink, IntraLink, LoopLink
- Filters: MinLen=4

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science



The raw data was peak picked in Mascot Distiller. Crosslinked precursors often have charge of 3+ or more, which means the product ions could also be highly charged. It's best to decharge the fragments by choosing to output fragment masses as MH+.

Setting up the crosslinking method is straightforward. DSS has a Unimod definition titled Xlink:DSS, and we only want the (K) specificity for lysine-lysine links. The authors determined only one of the three monolinks is possible, in this case the water quenched 156 Da monolink. This has the code W in the Unimod definition.

The accessions in the crosslinking method act as selectors and instruct Mascot to choose only the listed proteins for crosslinking. The sequences are in SwissProt, HOP2_ARATH and MND1_ARATH, so we can just list them here and select SwissProt as the search database when the search is submitted.

The crosslinked products to consider are protein interlinking and intralinking, and peptide looplinking. Finally, allow the beta peptide to be as short as 4 residues. If no MinLen is specified, Mascot filters out alpha and beta peptides shorter than MinPepLenInSearch, which is 7 by default.

Edit Crosslinking Method: PDX001538 Administrator Logout

Name
 Name: PDX001538 Description:

Method
 Strategy: Brute-force
 InterLink: ☒
 IntraLink: ☒
 LoopLink: ☒

Linkers
 Linker: Xlink:DSS (K)
 Monolink: ☒
 DoesNotPairWith: Xlink:DSS (K), Xlink:DSS (Protein N-term)

Accessions
 Database name: SwissProt
 Accession: MND1_ARATH
 Database name: SwissProt
 Accession: HOP2_ARATH

Filters
 Add parameter

Settings
 Add parameter

Save changes Cancel

Restrict pairing of linker specificities. The optional element can be used to restrict links between certain pairs of specificities. If there are no does_not_pair_with elements, a linker can connect any two linker specificities. Although it seems redundant, it is necessary to define both senses for each restriction. That is, if the linker element for K contains a does_not_pair_with element for Protein N-term, then the linker element for Protein N-term must contain a does_not_pair_with element for K.

Linker, linkable sites, monolinks

Protein selectors

MASCOT : Crosslinking

© 2020-2023 Matrix Science



The crosslinking method editor is shown here. Monolinks are toggled in the Monolink field. In the Accessions section you can optionally specify a database name and accession. If there is no database name, Mascot chooses proteins for crosslinking by accession, regardless of database. The filters and settings sections contain simple key-value parameters.

Interlinked HOP2-MND1 complex

The screenshot shows a web browser window with two tabs: 'Mascot search engine | Protein' and 'Mascot configuration'. The address bar shows 'localhost:8707/mascot/...'. The page title is 'Edit Crosslinking Method:PDX001538'. The user is logged in as 'Administrator' with a 'Logout' link. The form has a 'Name' field with 'PDX001538' and an empty 'Description' field. Below is a 'Method' section with a tab for 'XML'. The XML content is displayed in a text area, showing a configuration for a brute-force search strategy with specific linker and monolink settings. At the bottom, there are 'Update method', 'Save changes', and 'Cancel' buttons, along with a 'Help Window' placeholder.

Property	Value
xml	<pre><msm:method description="" name="PDX001538" strategy="Brute-force"> <msm:linkers> <msm:linker ModFileName="" link=DSS (K)> <msm:monolink>W</msm:monolink> </msm:linker> </msm:linkers> <msm:accessions> <msm:accession DatabaseName="SwissProt">PND1_AUTH</msm:accession> </msm:accession> </msm:accessions> </msm:method></pre>

MASCOT

: Crosslinking

© 2020-2023 Matrix Science

**MATRIX
SCIENCE**

The XML displays the raw method. It can be copied and pasted to other Mascot Server installations, shared with collaborators or Matrix Science support etc.

Interlinked HOP2-MND1 complex

- **Peak pick in Distiller (done)**
- **Set up a crosslinking method (done)**
- **Submit from Distiller**
 - Database: contaminants, SwissProt, *A. thaliana* taxonomy
 - Choose crosslinking method in search form
- **Open full report in browser**
 - Distiller support for displaying crosslinked matches will be added in a future release

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science



So, the raw data has been peak picked and we have a crosslinking method. The search is submitted as normal from Distiller. Since the two proteins are in SwissProt, I've chosen to search SwissProt with *A. thaliana* taxonomy for convenience. Also, remember to include a contaminants database. Choose the right crosslinking method in the search form.

Once the search finishes, choose to open the full report in browser. At the moment, Distiller 2.8 shows the crosslinked peptides in the tree. It can't show the peptide view for them though, so when you click on the match it'll open a browser windows with the peptide view from Mascot instead.

MASCOT Search Results

User :
 E-mail :
 Search title : 20140603_QEx3_RSLC4_Rampl_Mechtler_IMP_shotgun_DSS12.raw
 MS data file : 20140603_QEx3_RSLC4_Rampl_Mechtler_IMP_shotgun_DSS12.temp.mgf
 Databases : 1: contaminants 20090624 (262 sequences; 133,770 residues)
 2: SwissProt 2020_01 (561,911 sequences; 202,173,710 residues)
 Taxonomy : 1: (none)
 2: Arabidopsis thaliana (thale cress) (15,940 sequences)
 Timestamp : 11 May 2020 at 13:20:48 GMT

Re-search ☒ All ☐ Non-significant ☐ Unassigned [\[help\]](#) Export As XML

Search parameters

Score distribution

Modification statistics for all protein families

Modification	Delta	Type	Site	Total matches
Xlink:DSS[W]	156.078645	variable	K	465
Oxidation	15.994915	variable	M	340
Xlink:DSS	138.06808	crosslink	K - K	218
Xlink:DSS	138.06808	looplink	K - K	140
Carboxymethyl	58.005479	fixed	C	21

Legend

Protein Family Summary

Format Significance threshold p < 0.05 Max. number of families AUTO [\[help\]](#)
 Display non-sig. matches ☐ Min. number of sig. unique sequences 1
 Dendrograms cut at 0
 Preferred taxonomy All entries

Sensitivity

Tarnot

MASCOT : Crosslinking © 2020-2023 Matrix Science **MATRIX SCIENCE**

The Protein Family Summary report has been extended to display crosslinked search results. The modification statistics table gives counts of discovered intact links, monolinks and looplinks. Xlink:DSS[W] on the first row is the 156Da monolink enabled in the crosslinking method. The crosslink and looplink rows display the intact linker mass as well as the linked sites.

20140603_QEx3_RSLC4_Ramplic: X

localhost:8707/mascot/cgi/master_results_2.pl?file=.%2Fdata%2F20200511%2FF004610.dat

Format Significance threshold p < 0.05 Max. number of families AUTO [help]
 Display non-sig. matches ☐ Min. number of sig. unique sequences 1
 Dendrograms cut at 0
 Preferred taxonomy All entries

▼ Sensitivity

PSMs above homology 1180

Proteins (35) Report Builder Unassigned (9195) [s_permalink]

Protein families 1–10 (out of 35)

10 per page 1 2 3 4 Next Expand all Collapse all

Accession contains Find Clear

1	2::HOP2_ARATH	17682	Homologous-pairing protein 2 homolog OS=Arabidopsis thaliana OX=3702 GN=HOP2 PE=1 SV=1
2	2::MND1_ARATH	17505	Meiotic nuclear division protein 1 homolog OS=Arabidopsis thaliana OX=3702 GN=MND1 PE=1 SV=1
3	1::IPI:CON_Trypsin SWISS-PROT:P...	202	Trypsin - Sus scrofa (Pig).
4	1::IPI:CON_00009865.2 SWISS-PR...	75	Tax_Id=9606 Gene_Symbol=KRT10 Keratin, type I cytoskeletal 10
5	1::IPI:CON_00131368.3 SWISS-PR...	62	Tax_Id=10090 Gene_Symbol=Krt5a Keratin, type II cytoskeletal 6A
6	2::TBA2_ARATH	53	Tubulin alpha-2 chain OS=Arabidopsis thaliana OX=3702 GN=TUBA2 PE=2 SV=2
7	1::IPI:CON_00698900.1 SWISS-PR...	42	(Bos taurus) Actin, cytoplasmic 1
8	2::EDL18_ARATH	41	Sugar transporter ERD6-like 18 OS=Arabidopsis thaliana OX=3702 GN=SF22 PE=2 SV=1
9	1::IPI:CON_IPI00121462.1 SWISS-...	37	Tax_Id=10090 Gene_Symbol=Krt23 Keratin, type I cytoskeletal 23
10	1::IPI:CON_IPI00311493.5 SWISS-...	36	Tax_Id=10090 Gene_Symbol=Krt18 Keratin, type I cytoskeletal 18

10 per page 1 2 3 4 Next Expand all Collapse all

Mascot: <http://www.matrixscience.com/>

MASCOT

: Crosslinking

© 2020-2023 Matrix Science

**MATRIX
SCIENCE**

Protein inference is not affected by interlinking. Because HOP2_ARATH and MND1_ARATH don't share any significant peptide matches, protein inference puts them in different families, regardless of whether there are any interlinked peptide matches.

Notice also the small amounts of evidence for other Arabidopsis proteins in the sample, despite the high levels of purification.

-0.14	3	-	88	6.1e-09	►1	U	K.SLQSNLTLEEIQEKDAKLRK.E + K17<-Xlink:DSS[I]->K20	Looplinked peptide
-0.15	3	-	106	1.2e-10	►1	U	K.SLQSNLTLEEIQEKDAKLRK.E + Xlink:DSS[W]	Monolinks
0.39	3	-	123	3e-12	►1	U	K.SLQSNLTLEEIQEKDAKLRK.E + Xlink:DSS[W]	Intralinks in HOP2_ARATH
-7.75	3	-	49	7.5e-05	►1	U	K.EMEEKLVKLREGITLVRPEDK.K	
-2.16	1	1	59	3.2e-06	►1	U	K.TAVQKALDSLADAGK.I K5<-Xlink:DSS->K2 K.QKIYIAR.Q	
-2.09	1	1	34	0.00062	►1	U	K.TAVQKALDSLADAGK.I K5<-Xlink:DSS->K2 K.QKIYIAR.Q	
8.72	0	1	46	5.4e-05	►1*	[2:;MND1_ARATH]LES	DLQGSNK K10<-Xlink:DSS->K8 K.LQEQLQEKK.K	Interlink to MND1_ARATH

MASCOT

: Crosslinking

© 2020-2023 Matrix Science

MATRIX SCIENCE

Expand the second protein hit, HOP2_ARATH. The peptide match table displays crosslinking data in a compact form, and I’ve selected a few example rows here.

On the first line is a looplinked peptide. A looplink is an intact link between two residues in the same peptide. Here the intact link is between lysines K17 and K20. Looplinks use the symbol [I] in brackets to indicate it’s an intact link.

The next two rows show matches to the same peptide sequence, this time with the W monolink. Monolinks are variable modifications generated from the linker definition. With cleavable linkers, this is usually the only crosslinked product of interest, since monolinks represent the linker fragments. Monolinks can also be used for modelling dead-end or quenched links, such as the water quenched monolink here.

Further down, there are three matches with an intact crosslink. The first two are protein intralinks, which means it’s an intact link between two non-overlapping peptides within a single protein.

The last line shows a match to a protein interlink. An interlink is an intact link between two peptides from different proteins. The alpha peptide is in MND1_ARATH and the beta peptide in HOP2_ARATH.

The scores for all the matches are standard Mascot scores, and the identity and homology thresholds are calculated using the same statistical model as for linear peptide matches.

Validating intact crosslinked matches

- **Target-decoy searching?**

- Works with cleavable linker, because linker fragments are just variable mods

- **Intact linker**


- Sample has just two purified proteins
- Vast majority of peptide matches will cluster in the two proteins
- Too few sequences for meaningful FDR estimate

Which of the matches on the previous slide are correct? The usual approach is to do a target-decoy search, which allows you to estimate the false discovery rate. This approach works if you're using a cleavable linker. The linker fragment is just a variable modification and all the peptide matches are ordinary linear peptides., so standard FDR estimation techniques work well in this case.


However, when using intact linkers, the search space is often too small for a meaningful estimate of the FDR. In the present case, the sample is expected to contain only the two purified proteins. If you set up a two-sequence database and search ten thousand MS/MS spectra against it, no statistical model can give a meaningful estimate of the FDR. In this case, I've used as the target database the fifteen thousand *A. thaliana* proteins in SwissProt, plus a couple hundred contaminant sequences, which allows incorrect matches to distribute a bit more evenly. Even so, the peptide matches completely saturate the available peptide sequences from the two target proteins, and there are few matches in other proteins.

The Mascot statistical model is robust and will give you reasonable separation between correct and incorrect matches, even in these adverse conditions. Nonetheless, it's prudent to keep a sceptical mind. Let's look at a few example spectra.

All matches to this query

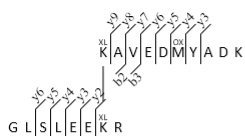

Score	Mr(calc)	Delta	Sequence
74.3	2253.1198	-0.0008	KAVEDMYADK K1<-Xlink:DSS->K7 GLSLEEK
22.1	2253.1198	-0.0008	KAVEDMYADK K1<-Xlink:DSS->K8 RGLSLEEK
1.5	2253.1198	-0.0008	KAVEDMYADK K10<-Xlink:DSS->K7 GLSLEEK

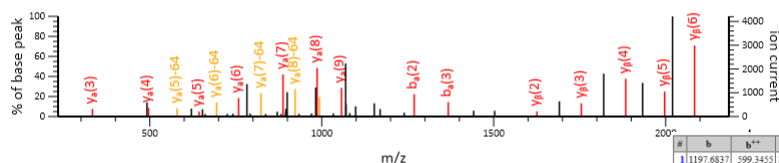

233.18 to 2184.01

Label all possible matches ☐ Label matches used for scoring ☒

Monoisotopic mass of neutral linked peptide Mr(calc) : 2253.1198
 Monoisotopic mass of neutral free alpha peptide Mr(calc) : 1184.5383
 Monoisotopic mass of neutral free beta peptide Mr(calc) : 930.5134
 Fixed modifications: Carbokymethyl (C) (apply to specified residues or termini only)
 Intact link : Alpha K1<-Xlink:DSS->K7 Beta
 Variable modifications:
 Alpha M6 : Oxidation (M), with neutral losses 0.0000 (shown in table), 63.9988
 Ions Score: 74 Expect: 5.5e-08
 Matches : 19/218 fragment ions using 36 most intense peaks ([help](#))

#	b	b ⁺⁺	b ⁺	b ⁺⁺⁺	b ⁰	b ⁰⁺⁺	Seq.	y	y ⁺⁺	y ⁺	y ⁺⁺⁺	y ⁰	y ⁰⁺⁺	#
1	1197.6837	599.3455	1180.6572	590.8322			K							10
2	1268.7089	634.8641	1251.6943	636.3508			A	1067.1507	529.2290	1040.4241	520.7157	1039.4401	520.2237	9
3	1367.7893	684.3983	1350.7627	675.8850			V	986.4135	493.7104	969.3870	485.1971	968.4030	484.7051	8
4	1496.8318	748.9196	1479.8053	740.4063	1478.8213	739.9143	D	887.3451	444.1762	870.3186	435.6629	869.3346	435.1709	7
5	1611.8588	806.4330	1594.8322	797.9198	1593.8482	797.4277	M	758.3025	379.6549	741.2760	371.1416	740.2920	370.6496	6
6	1758.8942	879.9507	1741.8676	871.4375	1740.8836	870.9454	D	643.2756	322.1414	626.2490	313.6282	625.2650	313.1362	5
7	1921.9575	961.4824	1904.9310	952.9691	1903.9470	952.4771	V	496.2402	248.6237	479.2136	240.1105	478.2296	239.6185	4
8	1992.9946	997.0010	1975.9681	988.4877	1974.9841	987.9957	A	333.1769	167.0921	316.1503	158.5788	315.1663	158.0868	3
9	2108.0216	1054.5144	2090.9950	1046.0011	2090.0110	1045.5091	D	262.1397	131.5735	245.1132	123.0602	244.1292	122.5682	2
10							K	147.1128	74.0600	130.0863	65.5468			1

#	b	b ⁺⁺	b ⁺	b ⁺⁺⁺	b ⁰	b ⁰⁺⁺	Seq.	y	y ⁺⁺	y ⁺	y ⁺⁺⁺	y ⁰	y ⁰⁺⁺	#
1	58.0287	29.5180					G							8
2	171.1128	86.0600					L	2197.1056	1099.0565	2180.0791	1090.5432	2179.0951	1090.0512	7
3	258.1448	129.5761			240.1343	120.5708	S	2084.0216	1042.5144	2066.9950	1034.0012	2066.0110	1033.5091	6
4	371.2289	186.1181			353.2183	177.1128	L	1996.9895	998.9984	1979.9630	990.4851	1978.9790	989.9931	5
5	500.2715	250.6394			482.2609	241.6341	E	1883.9051	942.4564	1866.8789	933.9431	1865.8949	933.4511	4
6	629.3141	315.1607			611.3035	306.1554	E	1754.8629	877.9351	1737.8363	869.4218	1736.8523	868.9298	3
7	2080.0154	1040.5114	2062.9889	1031.9981	2062.0049	1031.5061	K	1625.8263	813.4138	1608.7937	804.9005			2
8							R	175.1190	88.0631	158.0924	79.5498			1

MASCOT


: Crosslinking

© 2020-2023 Matrix Science

MATRIX
SCIENCE

Here's a Peptide View report of a crosslinked match with excellent fragmentation from both alpha and beta, with score 74. Further down the page, there is a fragmentation table for the alpha peptide and another for the beta peptide. The fragment masses have no overlap between the two, which shows both peptides have independent evidence.


At the bottom of the page, you can see the query also matches a slightly different beta peptide. However, the rank 2 score is much lower, only 22, and there's no ambiguity in the alpha peptide, so we can be confident the top match is the correct match.

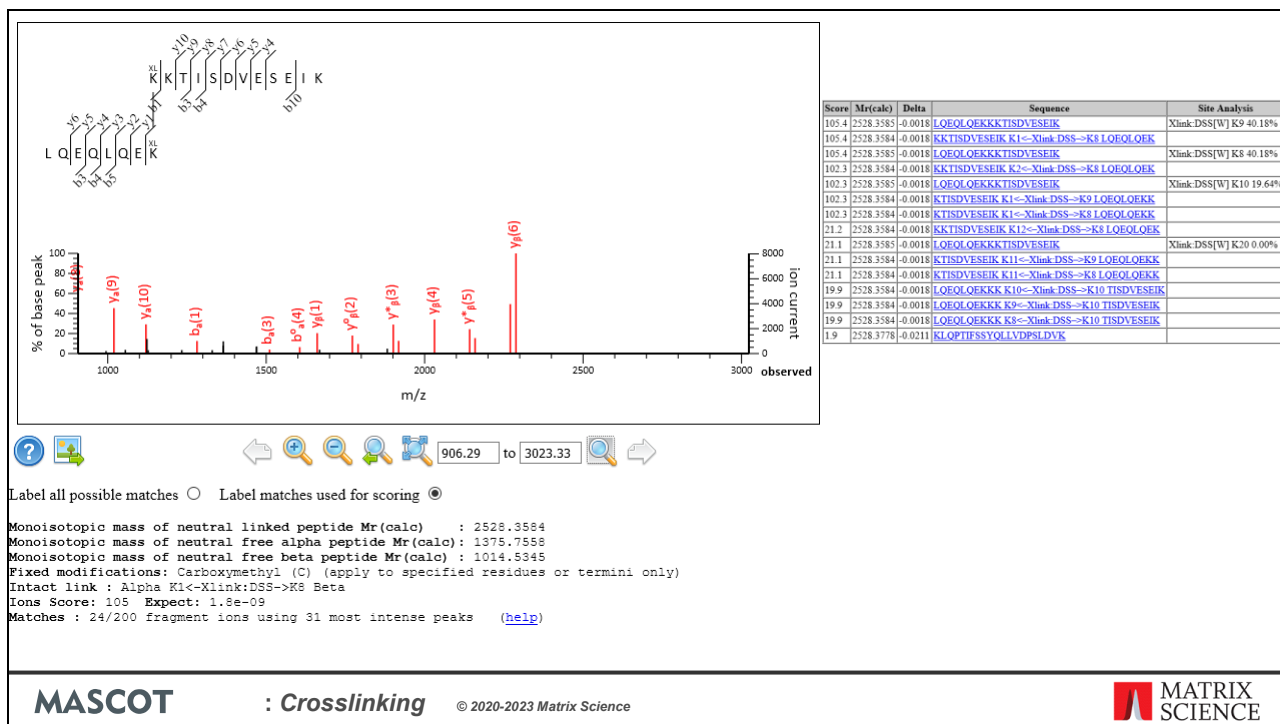

Here is another crosslinked peptide match with a high score and pretty nice fragmentation. At such a high score, 105, it looks very promising, but I have to dash your hopes.

Here are the matches at the other ranks. It turns out, beta and alpha are sequential in the protein sequence. The concatenated beta-alpha sequence with a water quenched monolink has exactly the same score and equally convincing fragment matches. The location of the potential monolink or crosslink is also in question. The MS/MS spectrum has no evidence to differentiate between the alternatives, so choosing the correct match requires consulting structural data.

This also shows the importance of presenting an integrated report. If the search engine matched only crosslinked or only linear peptides, the ambiguity could be easily missed.

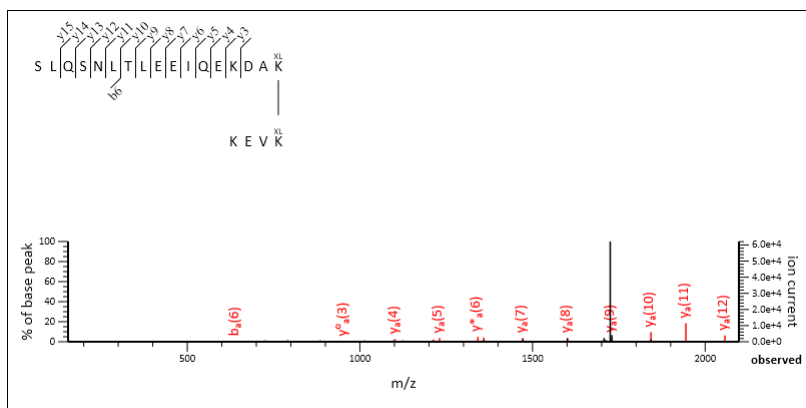

Score	Mr(calcd)	Delta	Sequence
87.2	2585.3799	0.0000	SLQSNLTLEEIQEKDAK K17<-Xlink:DSS->K4 KEVK
87.2	2585.3799	0.0000	SLQSNLTLEEIQEKDAK K17<-Xlink:DSS->K1 KEVK
84.4	2585.3799	0.0000	SLQSNLTLEEIQEKDAK K14<-Xlink:DSS->K4 KEVK
84.4	2585.3799	0.0000	SLQSNLTLEEIQEKDAK K14<-Xlink:DSS->K1 KEVK
84.4	2585.3799	0.0000	SLQSNLTLEEIQEK K14<-Xlink:DSS->K8 GVISQSVK


MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science

MATRIX
SCIENCE

Here's a crosslinked match where the identity of the beta peptide should be questioned. The alpha peptide has a nice sequence ladder, but there's no fragmentation from the beta. There's an almost as good a match at rank 5, where the beta peptide is completely different. The alternatives score almost the same because there are no alpha fragments that span the linker site. The two alpha peptides match exactly the same observed fragments.

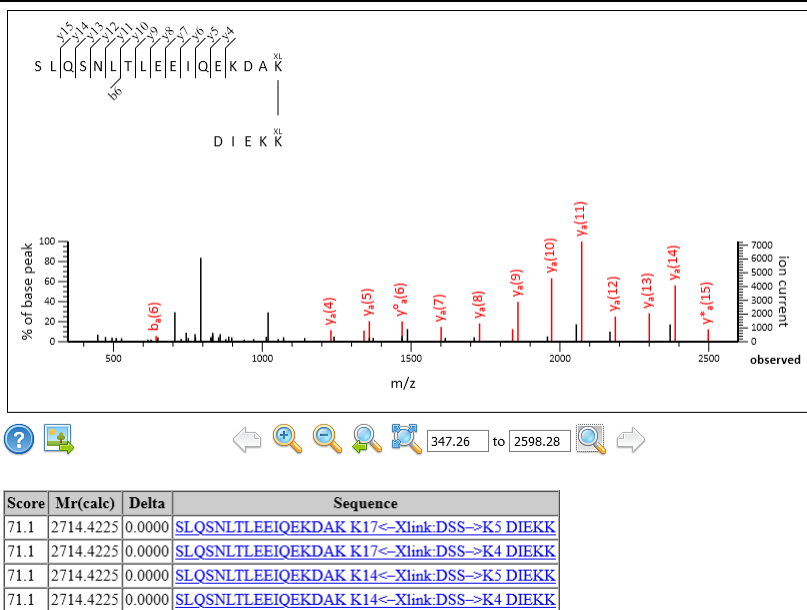

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science

**MATRIX
SCIENCE**

Finally, this crosslinked match is very similar to the previous one with no fragmentation from the beta peptide. But in this case, the search space has no other alternative for beta. If we assume the search space is complete, there's no reason to doubt the identity of the beta peptide, even if there are no beta fragments in the spectrum.

Visualising in xiVIEW

- **xiVIEW is a search engine independent, web-based visualisation tool**
- **Maintained by the Rappsilber laboratory**
- **<https://xiview.org>**

Other common ways to validate intact crosslinked matches is by comparing with existing annotations, and comparing with structural data. If the proposed crosslink would have a length longer than the linker's spacer arms, it's either a false positive or the structural data is incorrect. Mascot doesn't currently make use of structural data, so the validation must be done with external tools.

One way to visualise the results is by viewing them in xiVIEW. This is a search engine independent tool maintained by the Rappsilber laboratory.

Visualising in xiVIEW

- **xiVIEW CSV format**

- Choose xiVIEW CSV format in Mascot export form
- Also export peak lists as MGF in Mascot query order
- Export protein sequences as FASTA
 - Not needed if accessions are in sp|Q9FX64|HOP2_ARATH format

- **Or mzIdentML format**

- Under Protein Information, check Protein Sequence
- No need to check Matched Fragment Ions
- Also export peak lists as MGF in original order

- **Upload the data to <https://xiview.org>**

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science




You can export the results in two different formats for xiVIEW, either the xiVIEW CSV or mzIdentML format. xiVIEW documentation recommends mzIdentML.

If you use the CSV format, export also the peak lists as MGF and the protein sequences as FASTA. Make sure you export MGF in Mascot query order so they match with the CSV file.

If you use the mzIdentML format, enable Protein Sequence under Protein Information. This ensures protein sequences are saved in the mzIdentML file. Also export also the peak lists as MGF. In this case, make sure you export them in original order so they match with the order in the mzIdentML file.

Upload the files to xiview.org.

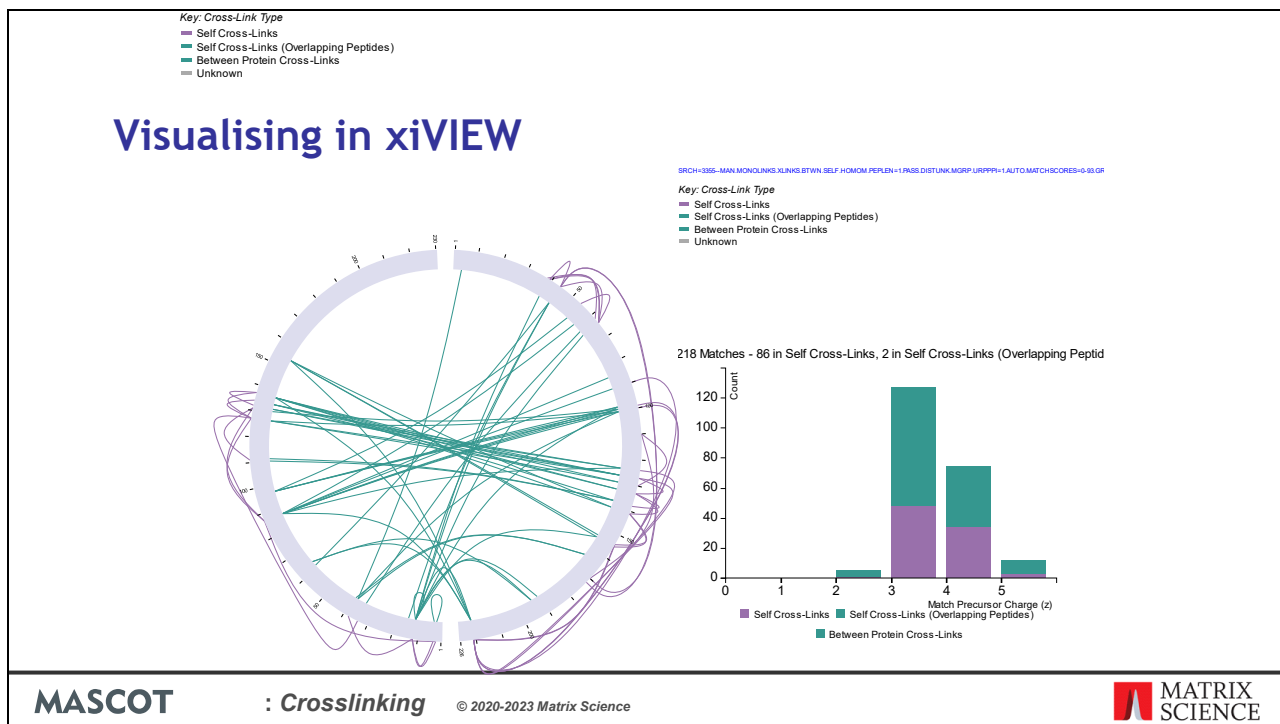

Here's an example circular plot of the crosslinks between the two proteins, based on the Mascot search data. xiVIEW can also plot histograms like this one showing the crosslinked match charge states. As expected, nearly all of the crosslinked matches have charge at least 3+ or higher.

Unfortunately, there is currently no crystallography data for the HOP2-MND1 complex from *Arabidopsis thaliana*, so I can't show you the 3D view.

Visualising in xiVIEW

	A	B	C	D	E	F	G	H	I	J	K	
1	Id	Protein1	PepPos1	PepSeq1	LinkPos1	Protein2	PepPos2	PepSeq2	LinkPos2	Score	CrossLinkerModMass	R
2	q1645_p1_1	MND1_ARATH	6	GLSLEEK	7	HOP2_ARATH	189	DVKELK	6	16.18	138.06808	
3	q2520_p1_1	MND1_ARATH	6	GLSLEEK	7	MND1_ARATH	35	M15.994915GPK	4	15.98	138.06808	
4	q3722_p1_1	MND1_ARATH	6	GLSLEEK	7	MND1_ARATH	2	SKKR	3	16.33	138.06808	
5	q3723_p1_1	MND1_ARATH	6	GLSLEEK	7	MND1_ARATH	2	SKKR	3	16.07	138.06808	
6	q4017_p1_1	MND1_ARATH	4	KRGLSLEEK156.078645R	1	MND1_ARATH	1	M15.994915SKK156.078645	3	24.69	138.06808	
7	q4974_p1_1	MND1_ARATH	120	TEALTQLK	8	MND1_ARATH	31	ELEK	4	33.83	138.06808	
8	q5157_p1_1	MND1_ARATH	6	GLSLEEK	7	HOP2_ARATH	39	KTAVQK	6	20.99	138.06808	
9	q5158_p1_1	MND1_ARATH	6	GLSLEEK	7	HOP2_ARATH	39	KTAVQK	6	28.67	138.06808	
10	q5284_p1_1	MND1_ARATH	6	GLSLEEK	7	HOP2_ARATH	221	KRPR	1	24.18	138.06808	
11	q5371_p1_1	MND1_ARATH	6	GLSLEEK	7	HOP2_ARATH	189	DVKELK	6	34.83	138.06808	
12	q6018_p1_1	MND1_ARATH	6	GLSLEEK	7	HOP2_ARATH	39	KTAVQK	6	33.64	138.06808	

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science



The xiVIEW CSV format is a simple tabular format, which you can load in Excel or LibreOffice. Modifications like oxidation are encoded as a delta mass within the peptide sequence. The CSV file can be edited before uploading to xiVIEW, for example if you want to remove a putative crosslink because fails a quality check.

General advice for successful analysis

- Make sure your peak picking program decharges the fragment ions to MH⁺ or outputs the fragment charge state
- Make sure the search space is complete - include the proteomes of the target species or expression system
- Only include the necessary protein sequences in crosslinking, no more
- Always look at the fragmentation and the alternatives other than the rank 1 match

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science



Crosslinked peptides are larger than typical tryptic peptides and have higher charge states. The resulting fragments also have higher charge states and need to be decharged or have the charge state reported.

Unless the sample is a synthetic sample where there is no background proteome, it's important to include the proteome databases for the target species or expression system in database selection.

Minimize the number of theoretical crosslinked peptides to be considered by only including the necessary protein sequences in crosslinking method.

Interpretation of the results is complicated, and you need to look at the fragmentation and the alternatives other than the rank 1 matches. There may be other interpretations of the data that have a similar scores that need to be accounted for.

More examples on our website

- **Reference documentation:**
<https://www.matrixscience.com/help/crosslink.html>
- **February 2020 blog article “Disulfide bond characterisation”**
- **Heterofunctional linker example:**
https://www.matrixscience.com/help/xl_examples.html
- **December 2022 blog article “Peak picking intact crosslink spectra with Mascot Distiller”**
- **March 2023 blog article “Analyzing the disulfide bond structure of the SARS-Cov-2 Spike protein”**

MASCOT

: *Crosslinking*

© 2020-2023 Matrix Science



If you'd like to see more examples of crosslinked searches, we have some on our website.